Image Merge in Elements

1. Make sure you have the channel tabs activated. Do this by clicking the Edit menu, then Options. Select the "Appearance" settings. Make sure the "Show Channel Tabs" box is checked.

General	Appearance	
Appearance Open Next Save Next Macro Measurement User rights	General Genera	
	Toolbars	
	Layouts ✓ Show Layout Tabs	
	Zooming mouse wheel - scrolls through Z stack for nd images containing Z - zoom in/out for other documents	
	CTRL + mouse wheel - zoom in/out for all documents	
	CTRL + mouse wheel - zoom in/out for all documents Keep text size while zooming	
	CTRL + mouse wheel - zoom in/out for all documents Image: Keep text size while zooming Image: Allow zoom factors lower than best fill Multiple Windows Image: Allow zoom factors lower than best fill	

2. open or acquire the first image. In the lower left corner of the image is the channel tab. Right click this tab. If you have not configured color configurations for your system (optical configurations) you will need to select properties, then change the channel name and color to suit the properties of that image.

3. Right-Click again and select "Copy Selected Channel".



4. Choose or acquire a second image. Again if needed on this new image set the channel properties (step 2). Place the cursor over the image and right-click. Select "Paste as New Channel".

1	Jones V	a french
	Show LUT Intensity	
Sec.	Hide Scrollbars	
1.	View Channel In Color	Ctrl+Alt+Shift+C
<	Paste As New Channel(s)	>
20	Paste As Brightfield	
Section Contraction	Calibrate Document	
28 / Uberellingt	Rescaling Method	•
Automated Mea	I <u>m</u> age Properties	I, Alt+Enter

5. You should now have a single image with two channel tabs. You can now save this image to produce a composite. You can also repeat these steps and add more channels to the current merge. There is no realistic limit on the number of channels one image can have.



Austin Blanco Technical Instruments <u>Austin@austinblanco.com</u>